



Wax-Off™

- General Information -

Wax-Off™ Paraffin Repellent prevents wax from sticking to equipment and countertops and has a pleasant smell. Use Wax-Off™ on microtomes, countertops and laboratory equipment. Wax-Off™ can also be used as a removing agent for solidified paraffin.

-Instructions for Use-

By using Wax-Off™, you save time in routine cleaning. Wax-Off™ replaces the need for harsh solvents, such as xylene and keeps equipment in pristine condition.

1. Spray a small amount on a paper towel or gauze pad
2. Apply a thin film to desired surfaces.

- Packaging -

Catalog#	Volume
WF1000	4oz Spray Bottle, CS/4
WF1030	30mL Spray Bottle, CS/12

- Precautions -

Wax-Off™ is safe for use on wood, metal, and plastic surfaces. Wax-Off™ dissolves paraffin, therefore it is not recommended for use on microtome blades, blade holders or base molds. Wax-Off™ has an odor similar to that of Spearmint. Wax-Off™ is a combustible liquid. Store locked up, and out of reach of children. May cause skin and eye irritation. See SDS for further safety precautions.

