



PathnSitu

Biotechnologies

Fast Freeze- Cryo Gel

Water Soluble OCT Compound
for Frozen Tissue Sectioning

- | | | |
|--------------------------|--------|--------|
| <input type="checkbox"/> | #FFG01 | Clear |
| <input type="checkbox"/> | #FFG02 | Blue |
| <input type="checkbox"/> | #FFG03 | Green |
| <input type="checkbox"/> | #FFG04 | Yellow |

118 mL. (4 Fl. Oz.)
For Laboratory Use Only
Made in USA

Date Received: _____ Date Opened: _____

Directions

Use Frozen Section Compound to secure and encapsulate tissue specimens for sectioning.

1. Form a flat base layer on surface of object holder, freeze.
2. Add additional Compound to cover base layer.
3. Place specimen on top of base layer.
4. Add more medium to encapsulate and cover specimen.
5. Use heat extractor to freeze and form flat sectioning surface.



WARNING

HAZARD STATEMENTS

Causes skin irritation. Causes eye irritation. May cause damage to organs (liver) through prolonged or repeated exposure.

PRECAUTIONARY STATEMENTS - PREVENTION

Wash body thoroughly after handling. Wear protective gloves. Do not breathe dust or vapors.

PRECAUTIONARY STATEMENTS - RESPONSE

IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing. If eye irritation persists: get medical advice/attention. Call a doctor if you feel unwell.

IF ON SKIN: Take off all contaminated clothing and wash it before reuse. Rinse skin with plenty of water. If skin irritation or rash occurs: get medical advice/attention. Specific treatment (see first aid section on this label).

IN CASE OF FIRE: Use CO₂ dry chemical, or foam for extinction.

STORAGE & DISPOSAL

Store locked up. Dispose of contents/container in accordance with local/national/international regulations.

NOTICE: Read Safety Data Sheet for additional information.

PathnSitu Biotechnologies

USA

538, Selby Ln
Livermore, CA 94588
Ph: +1 925-218-6939

INDIA

CDC Towers, 3rd Floor, B-Block
Plot#10/8, Nacharam IDA,
Road#5, Nacharam,
Hyderabad Pin 500 076
Ph: 040-2701-5533/5544

www.pathnsitu.com