



N-15™

ALDEHYDE NEUTRALIZER POWDER

- #N15010 10 BOTTLES/CASE
- #N15020 10 BAGS/CASE
- #N15030 BULK, 30 BAGS/CASE
- #N15032 BULK, 32 BOTTLES/CASE

For In Vitro Diagnostics Use



4300 Emperor Blvd - 400 Durham, NC 27703

Tel: 1-877-846-5393 Fax: 1-877-817-1716

www.cancerdiagnostics.com **Made in USA**

HK0209 N15 Bottle and Bag Primary Container AWW-1275 Rev.0 2020.04.14

Date Received:

Date Opened:

Directions: Empty the entire contents of bottle/bag into 1 gallon/4 liters of spent 10% formalin or up to 4% glutaraldehyde. Secure lid on container and agitate until powder is completely dissolved. Let solution stand for 20 minutes. You may use the aldehyde test kit to test for residual aldehyde. Dispose of waste to sanitary sewer system in accordance with local POTW requirements.



DANGER **HAZARD STATEMENTS**

Harmful if swallowed. Harmful in contact with skin. Causes serious eye damage. Suspect of causing genetic defects.

PRECAUTIONARY STATEMENTS - PREVENTION

Wash body thoroughly after handling. Do not eat, drink, or smoke when using this product. Wear protective gloves and clothing. Wear eye and face protection. Obtain special instructions before use. Do not handle until all safety precautions have been read and understood.

PRECAUTIONARY STATEMENTS - RESPONSE

IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing. Immediately call a doctor. If exposed or concerned: get medical attention.

IF ON SKIN: Wash with plenty of water. Call a doctor if you feel unwell. Specific treatment, see first aid section on the SDS. Take off immediately all contaminated clothing and wash it before reuse.

IF SWALLOWED: Call a doctor if you feel unwell. Rinse mouth.

STORAGE & DISPOSAL

Store at room temperature. Store locked up. Dispose of contents/container to an approved waste disposal plant.

NOTICE: Read Safety Data Sheet for additional information.

EMERGENCY CONTACT: Chemtrec: 1-800-424-9300