



# Cleanerol™ Spray

## - Intended Use -

Use Cleanerol™ Spray to remove paraffin wax, Ink, blood, stains, mold, tar, grease, crayon marks, oil, fresh paints, adhesives, asphalt, roof coatings, cosmoline petroleum-based products, graffiti, chewing gum, nail polish, glues and more.

## - Packaging -

Catalog#	Volume & Qty
CL2525	4oz CS/4

## - General Information -

A natural citrus solvent-based stain and tissue dye remover that is non-caustic, non-acid, and non-corrosive. Works great on microtomes, countertops, markers, and more. It's aggressive cleaning abilities come from D-Limonene, a by-product of citrus fruits. It can be used safely on a wide variety of surfaces. Cleanerol™ Spray provides outstanding penetration and dispersal and removal of grease, oil, crayon, marker, ink, tar, adhesives, and asphalt products. Water rinsable and 100% biodegradable.

## - General Directions -

Be sure to test in an inconspicuous area before using on glass, acrylic, some plastics and especially painted surfaces. Spray product full strength or dilute as necessary to a dry surface. Direction varies based on materials listed below.

- Hard Surfaces:** Allow Cleanerol™ Spray to remain on surface until spot or stain appears loosened, then rinse with water to remove residue.
- Fabrics:** Saturate fabric with Cleanerol™ Spray. Rub into stain, wash article of clothing in washing machine.
- Carpeting:** Apply Cleanerol™ Spray sparingly and blot stain lightly; repeat with clean cloth. Repeat as necessary, then rinse with water.
- Stubborn Stains:** Apply Cleanerol™ Spray and allow to remain for 10 minutes, then rinse with water.

