



TriCide10™

- Intended Use - KEEP OUT OF REACH OF CHILDREN

It is a violation of Federal law to use this product in a manner inconsistent with its labeling. Rinse hard, non-porous food contact surfaces with potable water after application of product. This product is not for use on medical device surfaces.

*Kills HIV, HBV and HCV on pre-cleaned hard, non-porous surfaces/objects previously Soiled with blood/body fluids.

- Active Ingredients -

Alkyl (50% C ₁₄ , 40% C ₁₂ , 10% C ₁₆)	
dimethyl benzyl ammonium chloride	0.034%
Octyl decyl dimethyl ammonium chloride	0.026%
Didecyl dimethyl ammonium chloride	0.013%
Diocetyl dimethyl ammonium chloride	0.013%
Other Ingredients	99.914%
TOTAL	100.000%

- General Instructions -

For use as a hospital disinfectant, virucide*, deodorizer, cleaner:

1. Pre-Clean heavily Soiled Areas.
2. Apply this product undiluted to disinfect hard, non-porous surfaces with a sponge, brush, cloth, hand pump trigger spray device, coarse trigger spray device. For spray applications, spray 6-8 inches from surface. **Do not breathe spray.**
3. Treated surfaces must remain wet for 10 minutes.
4. Wipe dry with a clean cloth or allow to air dry.
5. Change solution daily or when visibly dirty.

- Special Instructions -

For cleaning and decontamination against HIV-1, HBV and HCV on surfaces/objects soiled with blood/body fluids.

Personal Protection	When handling items soiled with blood or body fluids use disposable latex gloves, gowns, masks, and eye coverings.
Cleaning Procedure	Blood and other body fluids must be thoroughly cleaned from hard, non-porous surfaces and objects before application of this product.
Disposal of Infectious Materials	Blood and other body fluids, cleaning materials and clothing must be autoclaved and disposed of according to federal, state and local regulations for infectious waste disposal.
Contact Time	Allow hard, non-porous surface to remain wet for 2 minutes to kill HIV and for 10 minutes to kill all other viruses and bacteria listed on the product label.

- To Kill Fungi -

Pre-clean heavily soiled surfaces. Apply product to hard, non-porous surfaces. Allow surface to remain wet for 10 minutes. Wipe surfaces (and let air dry).

- Mold/Mildew Control -

Apply product to hard, non-porous surfaces which will effectively inhibit the growth of mold and mildew and their odors. Repeat treatment every seven days, or more often if new growth appears.

- Packaging -

Catalog#	Volume	Qty.
TC10-12	32 oz. Spray	CS/12

